


HOW TO USE SWABS FOR ENVIRONMENTAL TESTING

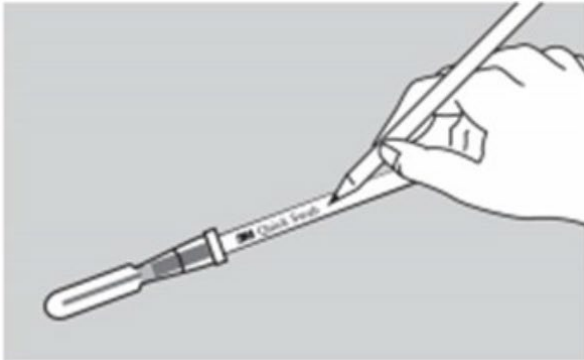


Clearly label swabs with a unique identifier.



Bend the top of the swab until you hear a snap.



Squeeze the bulb to dispense the broth onto the swab.

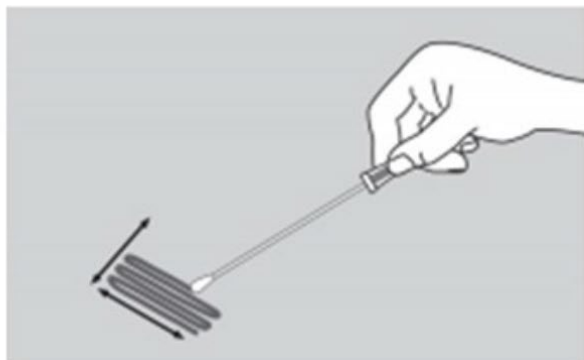

Sample an area no larger than 10 cm by 10 cm. Do not touch any other surfaces or hands with swab.

ATTENTION: Swabs must be refrigerated until used. They must be shipped cold, (2-8 degrees C) but not frozen in a cooler. They should be overnighted and received within 24 hours of testing to insure quality of results.

Please make sure that swabs are sampled and shipped prior to the expiration day on the package.

Please contact your account manager or call 402-334-7770 with questions

